

Pressing and Pelletizing




Fusion




XRF Accessories

## HANDBOOK OF SAMPLE PREPARATION & HANDLING

Pulverizing and Blending

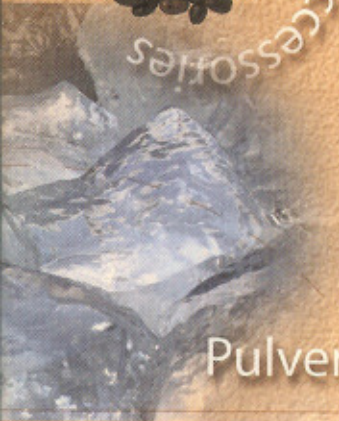

**SPEX**  
**CertiPrep**

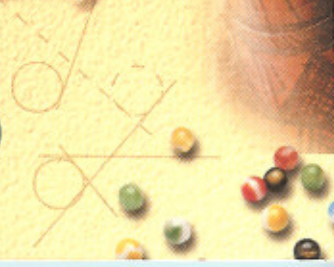

# **SPEX CertiPrep Handbook of Sample Preparation and Handling**

## **Sixth Edition Contributing Editors**

Ralph H. Obenauf, Ph.D  
Richard Bostwick  
William Fithian  
Mena DeStefano  
Joan D. McCormack  
David Selem

## **SPEX CertiPrep, Inc.**

203 Norcross Avenue  
Metuchen, NJ 08840

1-800-LAB-SPEX • 732-549-7144 • Fax 732-603-9647

e-mail: [sampleprep@spexcsp.com](mailto:sampleprep@spexcsp.com)

Website: <http://www.spexcsp.com>



# TABLE OF CONTENTS

Chemical Elements and Symbols .....	2
Periodic Chart of the Elements .....	3
How to Order .....	5
How to Use This Handbook .....	12

## SECTION I PRODUCT INFORMATION

### Chapter One Pulverizing and Blending

Introduction .....	14
Products for Pulverizing and Blending	
Boron Carbide Mortars and Pestles .....	15
Freezer/Mills .....	16
6750 Freezer/Mill .....	16
6800 Freezer/Mill .....	17
Freezer/Mill Accessories .....	17
6750 Accessories .....	18
6800 Accessories .....	19
Micro Hammer-Cutter Mill and Screens .....	21
Wig-L-Bug and Accessories .....	22
Mixer/Mills .....	24
5100 Mixer/Mill .....	24
Mixing and Grinding Containers for 5100 .....	25
8000D Mixer/Mill .....	28
8000 Mixer/Mill .....	30
Mixing and Grinding Containers (8000 Series) .....	31
Adapters and Smaller Vials (8000 Series) .....	36
Shatterboxes .....	39
8500 Shatterbox .....	39
8515 Shatterbox .....	40
8575 Flatterbox .....	41
Shatterbox Grinding Containers .....	42
Shatterbox Accessories .....	44
Nylon Sieves .....	45

### Chapter Two Pressing and Pelletizing

Introduction .....	48
Products for Pressing and Pelletizing	
3624B X-Press .....	49
3630 X-Press .....	50
3628 Bench Press .....	51
Manual Presses (SPEX CertiPrep-Carver) .....	52
Pellet Dies and Accessories .....	54
Sleeve-and-Plunger Sets for Pellet Dies .....	55
Replacement Pellets for Pellet Dies .....	56
Spec-Caps for Reinforcing XRF Pellets .....	57

### Chapter Three Fusion

Introduction .....	60
Products for Fusion	
SPEX CertiPrep-Claiss Automated Fusion Fluxers .....	61
SPEX CertiPrep-Claiss Fluxer Crucibles, Molds, and Accessories .....	64
Graphite Crucibles and Handling Equipment .....	67
Fusion Fluxes .....	70
Non-Wetting Agents .....	71

### Chapter Four XRF Accessories

Introduction .....	74
Products for XRF	
Spec-Caps for Reinforcing XRF Pellets .....	75
PrepAid Grinding Aid and KBr Pre-weighs .....	76
PrepAid Binders .....	77
X-Cells for XRF Liquid Samples .....	78
Thin Film for XRF Cell Windows .....	80
Certified Reference Materials (CRM Division) .....	83
Services Available .....	83
Products .....	84

## SECTION II TECHNICAL INFORMATION

### Chapter One

#### *Pulverizing and Blending*

Introduction .....	90
Grinding Container Materials .....	91
How to Select Grinding Containers .....	92
Charts	
Sample Container Materials Selection .....	93
Containers for Grinding and/or Blending .....	94
Key Points About Grinding .....	96
Summary of the Many Uses for Freezer/Mills .....	100
Products for Pulverizing and Blending	
Boron Carbide Mortars and Pestles .....	103
Comparison of Laboratory Mills—Chart C .....	104
Freezer/Mills .....	106
Guide to Freezer/Mill Capacity .....	108
Micro Hammer-Cutter Mill .....	111
Mixer/Mills .....	113
8000 Mixer/Mill Grinding Test Summary .....	118
Shatterboxes .....	119
Shatterbox Grinding Test Summary .....	123

### Chapter Two

#### *Pressing and Pelletizing*

Introduction .....	126
Preparing Powder Samples for XRF .....	127
Particle Size Effects .....	127
Pressing Sample Discs .....	129
Care and Maintenance of Pellet Dies .....	130
Set-up and Loading of Pellet Dies .....	131
Proper Use of Spec-Cap .....	132
Binders .....	133
Sleeve-and-Plunger Technique .....	135
Anti-sticking Agents .....	136
Pressing the Sample Disc .....	137
Sample Disc Removal .....	138
Use of Vacuum in Pellet Pressing .....	138
Sample Disc Handling .....	139
Trouble Shooting .....	140
Selecting a Pellet Die .....	141
Choosing a Laboratory Press .....	142
3630 Automated X-Press .....	142
3624B X-Press Motorized Hydraulic Press .....	143
3628 Bench Press .....	143
SPEX CertiPrep-Carver Manual Hydraulic Presses .....	144

### Chapter Three

#### *Fusion*

Introduction .....	146
When to Employ Borate Fusion .....	146
How to Prepare a Fused Sample .....	148
Selecting a Fusion Flux .....	149
Additives for Fusions .....	150
Products for Fusions	
Automated Fluxers .....	152
Graphite Crucibles .....	154

### Chapter Four

#### *XRF Accessories*

Introduction .....	156
Reinforcing XRF Pellets .....	157
X-Cells for XRF .....	157
Bubble-Free Cell Insert .....	158
Selecting Thin Film Windows for XRF .....	159
Thin Film Transmission Diagram .....	162
Chemical Resistance Tables .....	163
Bibliography .....	168
Trademark References .....	169

Catalog Number Index

Subject Index

Questionnaire for Test Grinding of Samples

Conversion Factors